



Fast Separation of Eleven Nitroaromatic Compounds on ZirChrom®-CARB

Clayton McNeff, Ph.D. and Kelly Johnson
ZirChrom Separations, Inc.

Technical Bulletin #280

This technical bulletin details the separation of eleven closely related nitroaromatic compounds, namely RDX, HMX, Nitrobenzene, 2-Nitrotoluene, Tetryl, 2,6-Dinitrotoluene, 4-Nitrotoluene, 1,3-Dinitrobenzene, 2,4-Dinitrotoluene, 2-amino 4,6-dinitrotoluene, 1,3,5-Trinitrobenzene. Our customers have reported that similar separations on silica-based phases produce run times as long as thirty minutes. Here we report a method on ZirChrom®-CARB at a column temperature of 125°C in under 4 minutes using a Metalox™ 200-C heater.

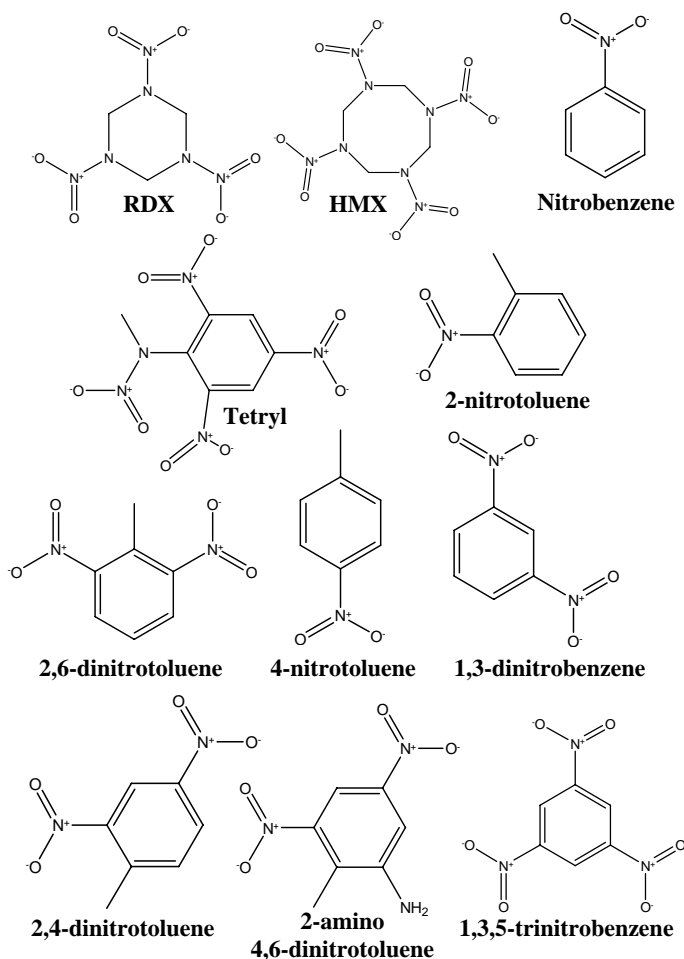


Figure 1: Structures of Explosives

Introduction

The rapid and accurate detection of nitroaromatic compounds is difficult due to the structural similarity of the compounds (see Figure 1). A new method developed at ZirChrom employs the unique temperature stability and surface chemistry of zirconia to achieve baseline resolution of these compounds in less than 4 minutes.

Experimental

A mixture of nitroaromatics was separated at 125°C using a ZirChrom®-CARB column (See Figure 2). The separation conditions were as follows:

Column: ZirChrom®-CARB,
150mm x 4.6mm i.d., (part#ZR01-1546)
Mobile phase: Isocratic Pre-mixed 35/15/50
Acetonitrile/Tetrahydrofuran/20mM ammonium carbonate pH 5.7, 10mM octylamine
Flow Rate: 2.0 ml/min.
Temperature: 125°C (Metalox™ 200-C Heater)
Detection: 254 nm
Inj. Volume: 1µl

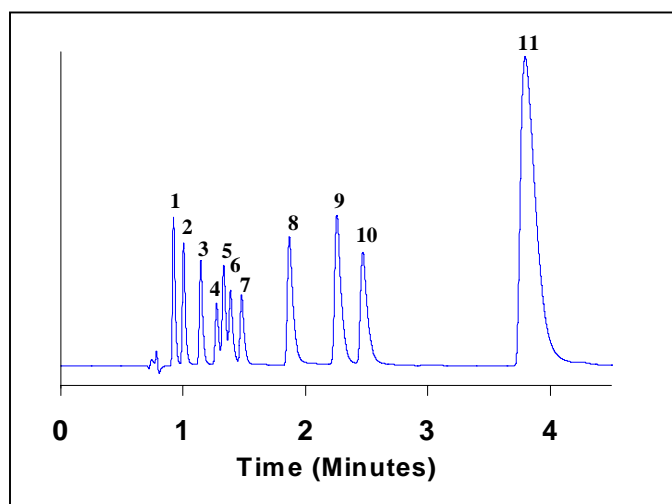


Figure 2: Separation of eleven nitroaromatics on ZirChrom®-CARB. 1=RDX, 2=HMX, 3=Nitrobenzene, 4=2-Nitrotoluene, 5=Tetryl, 6=2,6-Dinitrotoluene, 7=4-Nitrotoluene, 8=1,3-Dinitrobenzene, 9=2,4-Dinitrotoluene, 10=2-amino 4,6-dinitrotoluene, 11=1,3,5-Trinitrobenzene

This separation allows for clear identification and quantification of these compounds without the use of expensive MS detection. The separation is also completed using isocratic conditions, thus facilitating a more reproducible transfer from LC to LC.

Acknowledgments

We would like to thank Richard Burrows, Severn Trent Laboratories, for his many helpful discussions.

ZirChrom Separations, Inc.
617 Pierce Street, Anoka, MN 55303
1-866-STABLE-1
support@zirchrom.com

Visit www.zirchrom.com for more application notes using ultra-stable, high efficiency ZirChrom columns.